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
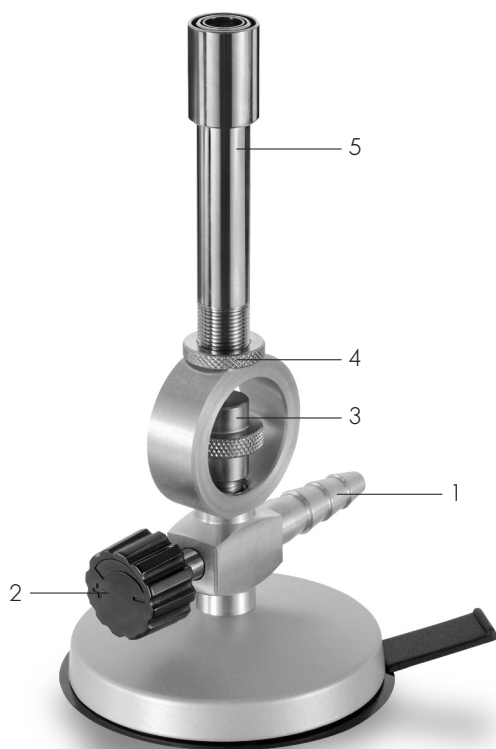
Funktion des Haffusses

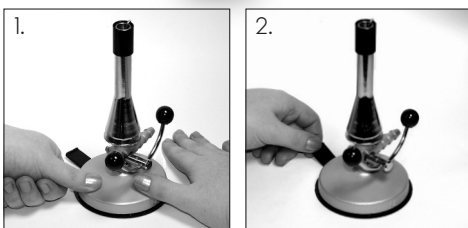Directions for use of Laboratory Burners

Validity:

The instructions are appropriate for:
Universal Burners.

Gas Pressure Range in Millibar

Nat. Gas: 18 to 25
L.P. Gas: 47,5 to 57,5 (use pressure regulator)
Town Gas: 7,5 to 15

Operating Instructions:

Fit gas tubing to inlet (1).
Use only DVGW approved laboratory tubing.
Secure with hose clip.
Open main gas tap.

– Turn needle valve (2) to left to open.
To select for required gas turn adjustable jet (3).

L.P. Gas: Open a little
Nat. Gas: Half open
Town Gas: Fully open

– Ignite.

Flame Type:

Loosen locking collar. (4)

Hard Flame: Screw burner tube (5) upwards.

Soft Flame: Screw burner tube (5) downwards.

After setting of tube secure with locking collar (4).

Safety Note:

Laboratory burners should only be used under constant supervision.
The main gas tap should be opened immediately before use and closed immediately after.
The Universal burner is supplied from the factory with secure needle valve, approved inlet and suitable sealers.

Function of safety base:

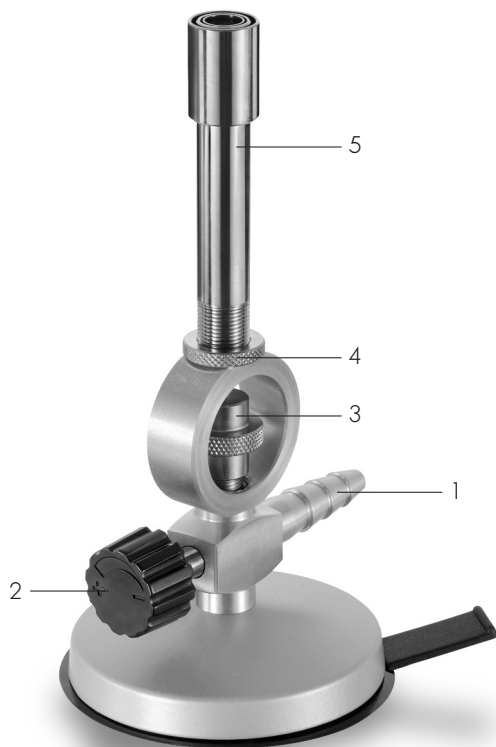
0. Screw safety base under laboratory burner base.

1. Press laboratory burner against clean and flat surface. (see picture)

2. To loosen draw tongue of safety base slightly upwards. (see picture)

If adhesion of safety base is reduced, clean with moist cloth.

Safety base available as spare part.

Function of safety base